Q800R3 DNA Shearing Sonicator





 Uses standard sample tubes

2 year warranty



APPLICATIONS:

- DNA shearing for Next-gen sequencing
- Chromatin shearing
- ChIP
- ChIP-seq
- RNA-seq
- Protein extraction
- Cell lysis

KEY INSTALLATIONS:

- Broad Institute
- NIH
- Harvard Medical School
- Broad Institute
- Memorial Sloan Kettering
- Max Planck Institute
- Ludwig Institute UCSD
- University of Cambridge
- Washington Univ. St. Louis
- California Institute of Technology
- Shanghai Inst. for Biological Sciences

Qsonica has over 45 years of experience manufacturing ultrasonic equipment with thousands of customers around the world. All products are designed and manufactured in the USA and include a full 2-year warranty.

The Q800R3 is our 2nd generation DNA and Chromatin shearing system. Improvements include an enhanced user-friendly design and quieter operation while remaining thousands of dollars less than the competition.

Samples can be sheared to a range of fragment sizes (150bp – 3kb). The system is compatible with commercially available sample tubes and is capable of processing up to 18 samples at one time. Multiple tube rack options can accommodate 50ul - 1ml sample volumes using a variety of standard sample tubes.

The Q800R3 is a complete package including a chiller to control temperature, digital operating system with programmable memory, high intensity ultrasonics and sound reducing enclosure.

Chromatin Shearing

Example protocols and results are based on customer feedback.

Sample Protocols:

Mammalian Chromatin

Cell Type: HEK 293T (2) 10cm dishes

70-80% confluent

Total Sample Volume: 300ul

Fixation Time: 1% Formaldehyde, 13 min

Sonicator Amplitude Setting: 70%

Sonication Pulse Rate: 15 seconds On,

45 seconds Off

Total Sonication On Time: 30 min

Yeast Chromatin

Cell Type: Wild type *S. pombe* cells grown to an OD of 1.3-1.5 in YEA.

Total Sample Volume: 300ul

Fixation Time: 1% Formaldehyde, 15 min

Sonicator Amplitude Setting: 100%

Sonication Pulse Rate: 20 seconds On,

40 seconds Off

Total Sonication On Time: 30 min

C. elegans Chromatin

Cell Type: *C. elegans* culture; Nuclear extract 1mg/ml

Total Sample Volume: 500ul

Fixation Time: 2% Formaldehyde, 20 min


Sonicator Amplitude Setting: 70%

Sonication Pulse Rate: 30 seconds On,

30 seconds Off

Total Sonication On Time: 20 min

Sample Results:

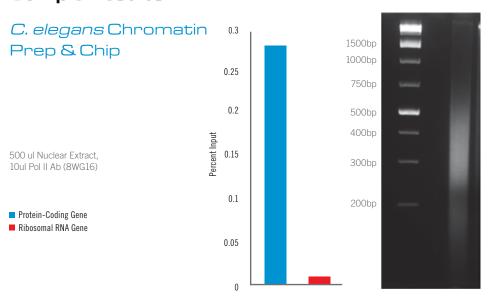

Mammalian Chromatin Replicates 2 Prep & Chip Chromosome 10* 6 Percent Input 2 300 bp H3K9me3 H3K9Ac H3K36me3 Н3 Chr10 Primer sequences:Forward - TCCTTCTCCCAACAATCAGCReverse - GATGTCGCTCCGAATCTTG Anithodies Used: H3K9me3 (abcamab8898), H3K9Ac (upstate 07-352), H3K36me3(abcam ah9050) H3 (ahccamah1791) * Average of two independent replicate chromatins in pane

Yeast Chromatin Prep & Chip



Sample protocols and publications for additional species and cell lines are available in the literature section of www.sonicator.com

Sample Protocols:

Bacterial Genomic DNA

Cell Type / Concentration: E.coli / 250ng

Sonication Pulse Rate: 15 seconds On, 15 seconds Off

Total Sample Volume: 200ul in 0.5ml tubes **Total Sonication On Time:** As indicated below

Sonicator Amplitude Setting: 20%

Human Skeletal Myoblast Genomic DNA


Cell Type / Concentration: LHCN-M2 Skeletal Myoblast / 200ng

Sonication Pulse Rate: 15 seconds On, 15 seconds Off


Total Sample Volume: 200ul in 0.5ml tubes **Total Sonication On Time:** As indicated below

Sonicator Amplitude Setting: 20%


Fragmentation Over Time

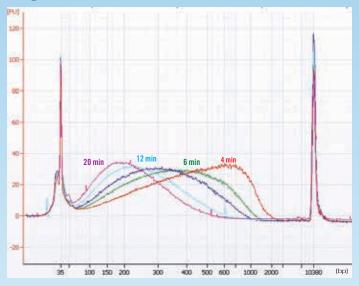


Fragmentation Over Time

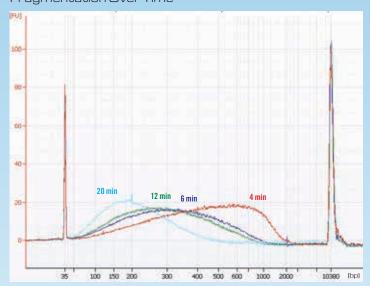

Testing Independent Replicates

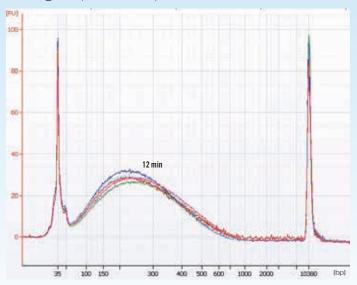

Fragmentation Over Time

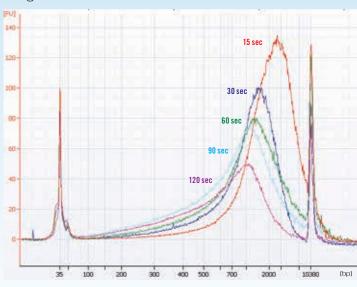



Sample Tube Rack Options for the Q800R3 System

#4256

8 tube holder (1.5mL Polystyrene tubes)



#425512 tube holder

12 tube holder (0.5mL PCR tubes)



#4262

18 tube holder (0.3mL PCR tubes)



Technical Specifications:

Power Rating: 750 Watts **Frequency:** 20 kHz

Programmable Timer: 1 second to 10 hours **Voltage:** 110V*, 50/60Hz

*Specify desired voltage for export

Dimensions:

Generator 8.00" W x 15.25" L x 8.50" H Enclosure 11.50" W x 12.00" L x 20.00" H Chiller 11.00" W x 13.00" L x 13.00" H

Additional information on our website:

- Sample Protocols
- Optimization tips
- Video
- Customer Feedback
- Recent Publications



Qsonica, LLC. 53 Church Hill Road, Newtown, CT 06470 Phone: 203.426.0101

E-mail: Info@Sonicator.com

